

Advanced Microscopy Services



Advanced microscopy techniques such as SEM (Scanning Electron Microscopy) (<https://www.eag.com/scanning-electron-microscopy-sem/>), TEM (Transmission Electron Microscopy) (<https://www.eag.com/tem-stem/>) and Dual Beam SEM (<https://www.eag.com/techniques/imaging/dual-beam-fib/>) are essential techniques to investigate sample microstructure, morphology, particle size, particle coatings and defects. These techniques often employ elemental mapping capabilities such as EELS (Electron Energy Loss Spectroscopy) (<https://www.eag.com/electron-energy-loss-spectroscopy-eels/>) and EDS (Energy Dispersive X-Ray Spectroscopy) (<https://www.eag.com/energy-dispersive-x-ray-spectroscopy-eds/>) which provide valuable information about elemental composition and location/distribution.

EAG offers a broad range and large installed base of different microscopy (<https://en.wikipedia.org/wiki/Microscopy>) tools and services to match your application, ranging from process development to failure analysis. In addition to providing high resolution imaging, our analytical capability makes us a unique partner that can help you during research, development, and analysis of failures.

Microscopy techniques can be used to characterize many types of defects. These include:


- Metal migration: FIB cross-section and investigation by STEM/EDS
- Voids: FIB cross-section and investigation by SEM
- Particles: Surface imaging or FIB cross-section to investigate size, chemistry and position in the layer stack
- Cracks and Delaminations: Small area FIB cross-sections or large area argon ion milled cross sections determine the location and of the delaminated interface. If needed, the interface chemistry can then be determined by TEM micro-analysis.
- Thickness and Uniformity: FIB cross-section and investigation by SEM

The types of materials tested using advanced microscopy techniques:

- Nanoparticles
- Alloys and metals
- Thin films
- Coatings on glass, silicon or carbon-based substrates
- Ceramics
- Composite materials
- IC devices
- Aluminum anodization


Industrial Applications of Microscopy

There has been a rapid emergence of new technologies which has enabled new industries such as facial recognition, autonomous driving, virtual reality, and 5G communication. These growth areas typically include FinFET, VCSEL, and III/V compound semiconductors. For our Advanced Microscopy team the materials indicate hetero-epitaxy, cubic and hexagonal lattices, where failures in devices often have complicated 3D structures.

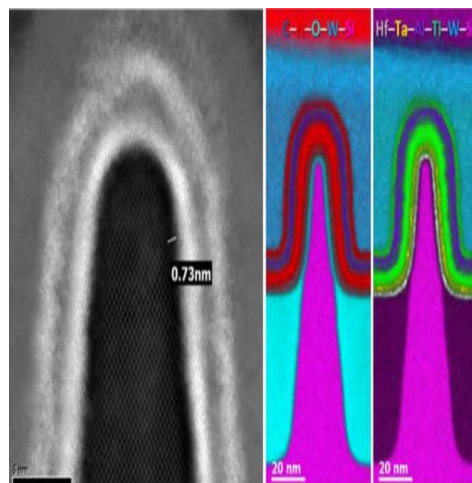


Recent development of FIB technology enables high-precision TEM sample preparation of 3D logic and memory devices. This example shows a HAADF image of a gate cut from a 22nm generation 3D FinFET device with $L_{\text{Gate}} \sim 30\text{nm}$ as seen in a recent paper by EAG Laboratories Advanced Microscopy team that was presented at Microscopy and Microanalysis 25 (S2): 690-691 (2019) (<https://www.cambridge.org/core/journals/microscopy-and-microanalysis/article/industrial-applications-of-electron-microscopy-a-shared-laboratory-perspective/B0FC0E2BCDC49A3ABF2601B66D96309D>) "Industrial Applications of Electron Microscopy: A Shared Laboratory Perspective"

Another interesting example is the picture below that shows a Aluminum Nitride/Silicon Carbide interface observed using HR-STEM imaging.

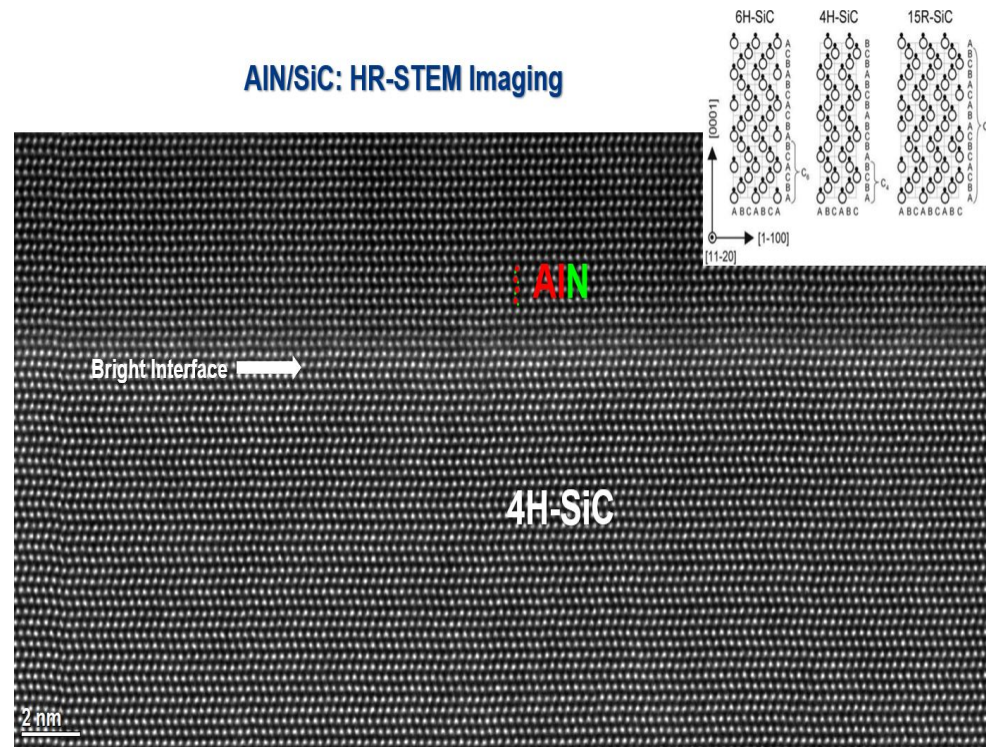

In the example below, we observe GaN Dislocation Typing: Burgers Vector Analysis with Large-Angle Convergent-Beam Electron Diffraction (LACBED)

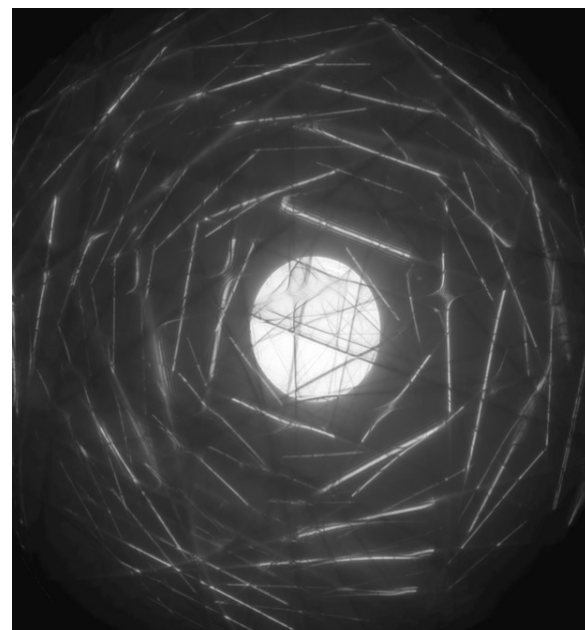

$$\begin{aligned}
 1u + 1v - 2t + 8w &= +9 \\
 0u - 1v + 1t - 4w &= -4 \\
 -2u + 4v - 2t + 6w &= +4 \\
 b &= [uvw], t = -(u+v) \\
 \text{Solution: } u &= \frac{2}{3}, v = -\frac{1}{3}, t = -\frac{1}{3}, w = 1 \\
 \text{Dislocation Burgers Vector} \\
 \mathbf{b} &= \frac{1}{3}[2\bar{1}\bar{1}3] \\
 &= \frac{1}{3}[2\bar{1}\bar{1}0] + \frac{1}{3}[0003] \\
 \mathbf{a+c} &\text{ mixed dislocation} \\
 \text{To verify the answer:} \\
 -1u + 0v + 1t - 5w &= -6 \\
 0u + 2v - 2t + 7w &= +7
 \end{aligned}$$

Would you like to learn more about our Advanced Microscopy services?

Please complete the form below to have an EAG expert contact you.

First Name*

Last Name*

Business Email*

Company Name*

Phone Number

Country
Please Select (reqd for ITAR compliance) ▾

How can we help you?*

How would you like us to contact you?*

- Phone
- Email
- No Preference

From time to time, we would like to share scientific content or EAG news that may be of interest to you. If you consent to our contacting you for this purpose, please check below.

I agree to receive communications from EAG Laboratories.*

You can unsubscribe from EAG communications at any time. To learn more, please see our Privacy Policy (<https://www.eag.com/privacy-policy/>).

By clicking below, you consent to allow EAG to store and process the personal information you submit.

SUBMIT

