



PROVIDING RESEARCHERS THE BEST INSTRUMENTATION
AT AN ECONOMICAL PRICE

- HOME
- PRODUCTS
- VIDEO TUTORIAL
- LEARN & EXPLORE
- TESTIMONIAL
- CUSTOMER PUBLICATIONS
- WHERE TO BUY
- CONTACT

[Products](#) > [IP Systems](#) > Basic IP System

InCyt Basic I/P Micro-Photometry System

- Ratiometric(Fura-2) imaging as fast as 1.5 ratio/sec. Single wavelength (Fluo-3) as fast as 15 pts/sec. Ratiometric photometry as fast as 6 ratio/sec. Single wavelength photometry as fast as 100 pts/sec.
- Ca++, pH, NO, Na.
- Other applications: FRET, GFP, vessel diameter, cell length, uncaging.
- Complete Turnkey systems.


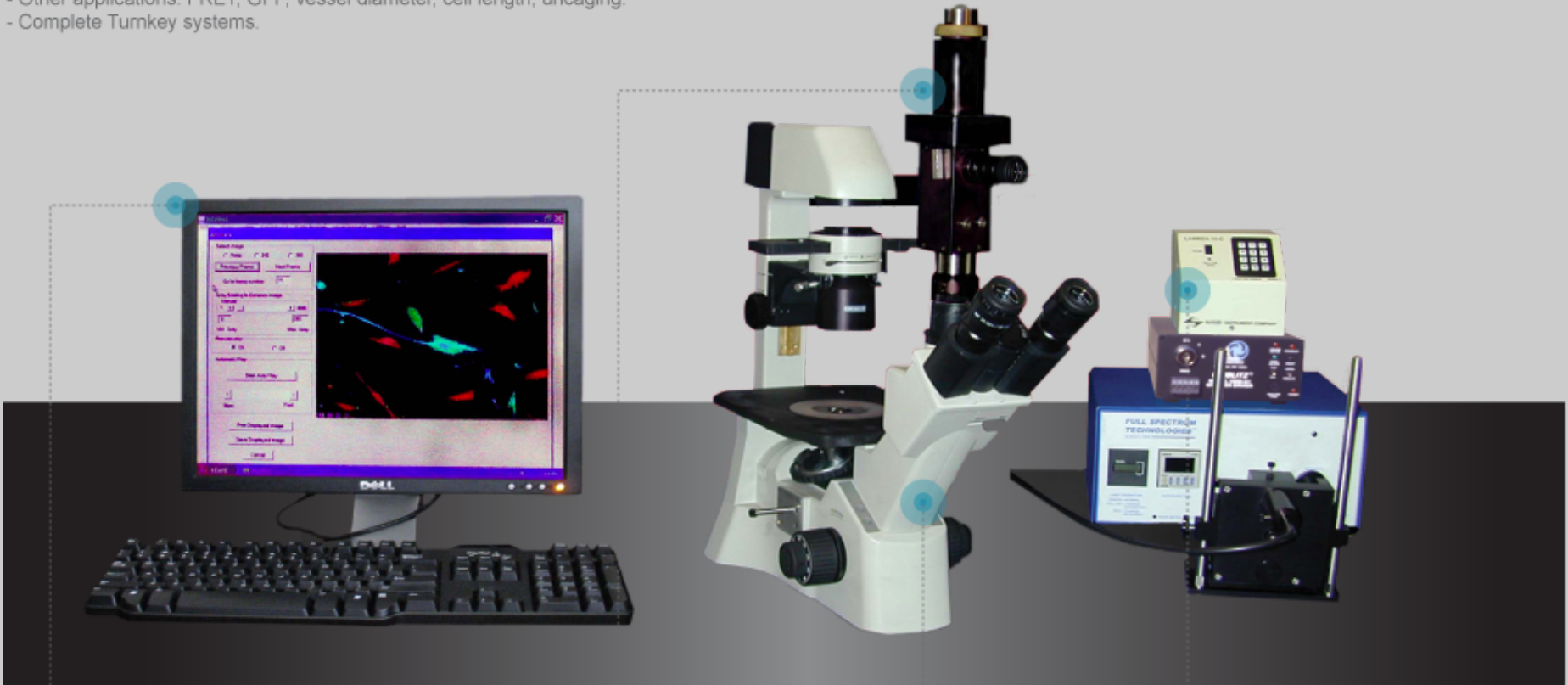


Image Acquisition & Analysis Workstation

State of art PC with Windows based proprietary InCyt software. Intuitive user interface guides the investigator through the experiment.

Photometer

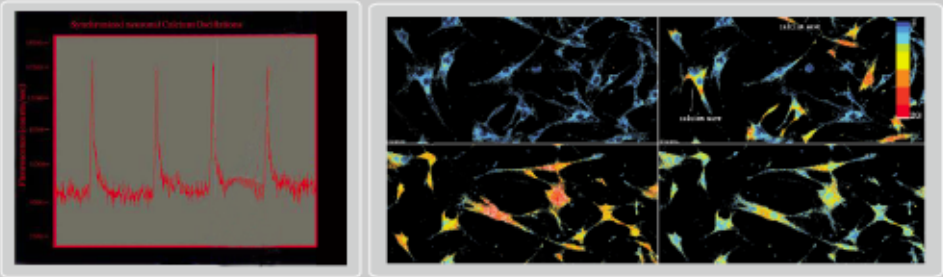

Exclusive design includes the following features:

- Photon counting for extraordinary sensitivity.
- Capture cell responses as fast as 10 milli-seconds.
- Variable aperture to isolate any area of the field of view.
- Computer-controlled shutter for PMT overload protection(optional).

Camera



Integrating 12 bit CCD video camera for low light level and low noise imaging.

Microscope

Accuscope 3032 Inverted Fluorescence Microscope with proprietary Groony fluorescence module.

Illumination System

Computer-controlled Sutter Lambda 10-b filter wheel with 120 watt Mercury-Xenon light source.

Complete Turnkey System

The InCyt Basic I/P™ Imaging System includes everything needed for calcium and other single excitation and ratiometric fluorophore experiments. The system includes: an inverted epi-fluorescence/phase contrast microscope, low-light level integrating CCD camera, computer controlled filter changer, Xenon UV/VIS arc lamp, image processing computer and data acquisition/analysis software.

Ease-of-use

The InCyt Basic I/P™ Imaging System was designed in a biomedical research laboratory by scientists for scientists. Menu selection appears in the sequence in which experiments are performed. Esoteric options that add complexity and are rarely used have been eliminated from the menus. The software follows standard Windows graphical-user-interface-protocol. The hardware has been tested for reliability and ruggedness, as well as, simplicity of set up and operation. New users can learn to use the system and be doing their first experiment in a few hours.

Data Collection, Analysis & Presentation

Data can be measured continuously on as many as 50 pre-selected regions of interest. The results can displayed in real time or images can be saved

for *post hoc* analysis. Saved images can be played back as an animated sequence—ideal for analysis of sub-cellular responses, intercellular communication and heterogeneity among the individual cells in a population. Ion concentration within cells can be converted to color using standard or custom pseudo-color tables. Image sequences can be presented as a montage with user-defined annotation for presentation and publication. Data generated by **InCyt Basic I/P™** are stored in a standard TIFF format and data files are stored as ASCII text.

Specifications

Camera Module	
Camera	Basler Scout Monochrome Integrating 12-bit CCD
Excitation Light Source	Sutter Instruments LB-LS17 light source with 175W Xenon arc lamp for both UV and visible light excitation.
Filter Changer	Tofra Wheel Filter Changer
Image Resolution	640 x 480 pixels
Camera Module	
PMT	Hamamatsu HC135-11 photon counter
Linearity (precision)	+/-1% from 0 to 20,000,000 counts/second
Maximum Dark Count (noise)	150 counts/second
Equivalent Noise Input (noise)	3 x 10 ⁻¹⁷ watts per second (measured at 400nm)
Spectral Sensitivity	300-650nm
Discriminator & Microprocessor	Built into PMT
Overload Protection	Computer-controlled shutter (optional)
Excitation	Sutter Instruments LB-LS17 light source with 175W Xenon arc lamp for both UV and visible light excitation.
Filter Changer	Tofra Wheel Filter Changer
Data Acquisition and Data Analysis Workstation	
CPU	Intel® 2nd generation Core™ i3 Dual Core Processor

RAM	4GB Non-ECC dual-channel 1000MHz DDR3 SDRAM
Hard Drives	250GB and 1TB 3.5" SATA 6Gb/s with 8MB DataBurst Cache
DVD ROM	16X DVD +/- R/W
Operating System	Windows 7
Monitor	20" SVGA

Data Capture and Display	
Maximum Speed	1.5 ratiometric points per second 15 single-wavelength points per second
Object Definition	Specify up to 50 user-defined regions for separate analysis
Data Collection Options	(1) Graph ion concentration in each cell during experiment. (2) Save images for later analysis and animated playback
Image Forma	TIFF
Data Storage Format	Tab-delimited ASCII

Other	
Calibration	From standard solutions or formula
Ions	Works with hundreds of fluorescent dyes for measuring: Ca ²⁺ , pH, Na ⁺ , K ⁺ , Mg ⁺² , H ₂ O ₂ , cyclic ADP Ribose, etc.

Data Presentation

- Animated playback of images
- Montages of selected images with user-defined annotation
- Graphs of ion kinetics in user-defined regions
- [For More Information...](#)

Corporate Headquarters: Dave Will
Service Department: Jess Luna
Research & Development: Eric Gruenstein
Marketing: Eric Vanderhoef